Application Note High-current Pulses for Battery Research



Introduction

There are many different ways of testing batteries and battery materials to determine their characteristics. Traditional methods include long-term cycling to determine cycle life and capacity fade, electrochemical impedance spectroscopy (EIS) to internal resistance, capacitance and other characteristics, potentiostatic (and galvanostatic) intermittent titration techniques (PITT/GITT) for diffusion rates, leakage rates, and self-discharge. There are many others but these form the basis for a large number of tests.

Recently, there has been an interest in rapid, higher current pulsing of batteries. This is different than what you get when you test using an Urban Drive Cycle where you have a series of pulses to simulate start/stop driving during a normal car ride. These pulses are on the order of 100 μ s to 100s of ms and can be used to stress a battery in ways that normal DC charging/discharging can't.

Gamry Instruments equipment has long been able to do these types of rapid, highcurrent pulses. One such example is the Reference 3000 potentiostat/galvanostat and Reference 30k Booster. Together, these two pieces of equipment let you pulse up to 30A on the order of 10s of microseconds. Now, just because our equipment can handle these types of demands doesn't mean your cell can. This application note is meant to show that our equipment can handle these demanding applications. It is up to you then to take it one step further and see if your cell can respond to these types of pulses.





Fig. 1. Reference[™] 3000 (left) with Reference[™] 30k Booster (right), with Test and Cal cell (green).

As a cell, we used our 30k Test and Cal cell. The 30k Test and Cal cell printed circuit board was carefully designed for:

- True four-terminal sensing on both the Test and Calibration sides of the Cell
- Minimal inductance in the current-carrying connections
- Minimal mutual inductance between the current-carrying and sensing circuits.

Gamry Instruments uses the Calibration side of the cell to galvanostatically measure impedance spectra at frequencies up to 300 kHz with less than 2° of phase error.

This Application Note shows results of our own research into the characteristics of our equipment.

Experimental

The tests were run on a Reference 3000 and Reference 30k Booster. A Gamry Reference 30k Test and Cal cell was connected to the booster in all tests. The cell connections used the standard cables supplied with the system: a 1 m currentcarrying cable and a 1.5 m sense cable.

Most of the testing was done on the calibration side of the Test and Cal cell. This side is a 200 m Ω resistor with a 2.5 A fast-blow fuse in the counter electrode lead. The fuse protects the 1 W resistor from damage caused by excessive current flow. The resistance of a sample fuse was measured to be 34 m Ω . Fast current pulses larger than the 2.5 A fuse rating will not blow the fuse if the pulses are short enough.



The Test side of the Test and Cal Cell is a 3 m Ω four-terminal resistor. It does not require a fuse because it can safely pass 30 A currents.

Many tests were performed at different CA (control amplifier) speed settings. The faster the CA speed is, the better time-resolution is achieved, but at the cost of stability in the amplifier. For a review of potentiostat design, see our Fundamentals of a Potentiostat Application Note.

Some of the tests required the additional use of a Tektronix 2024C digital oscilloscope in order to show CA response. This was achieved using the by attaching the oscilloscope to the E or I monitor BNCs on the rear of the potentiostat.

Results

2 A Pulses on 200 mΩ Cell, Using the Framework[™] Software

Fig. 2 below shows 2 A current steps lasting 200 µs each. Connections were made through the Reference 30k Booster, with the Reference 3000's 3 A full-scale range.

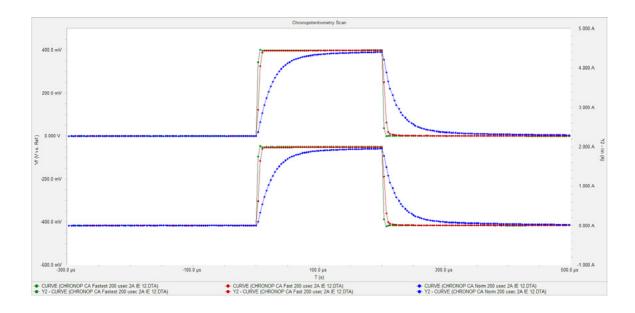


Fig. 2. The cell voltage is the upper trace, and cell current is the lower. Colors correspond to different control-amplifier speeds (CA speeds): green are CA speed **fastest**, (red) are CA speed **fast**, (blue) are CA speed **norm**.

Note the symmetry between the rising and falling waveforms. Charge lost from one edge of the pulse is recovered on the other edge.

4



Oscilloscope Data on 200 $m\Omega$ Taken from the E Monitor BNC

A BNC cable was connected between the Reference 3000 rear panel E Monitor output and the input of the oscilloscope.

Fig. 3 shows the voltage-versus-time curve of a 200 µs, 2 A pulse on the Calibrationside resistor. The graph was recorded in CA speed **fast**.


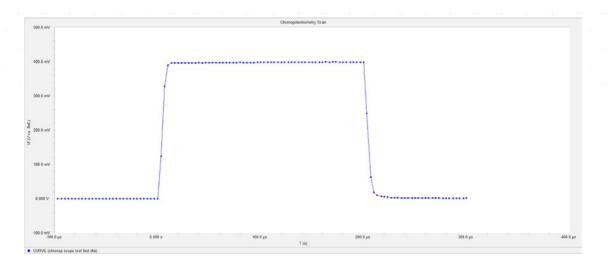


Fig. 3 Data recorded through the E Monitor connector.

The oscilloscope-screen capture (Fig. 4) shows a very similar waveform.





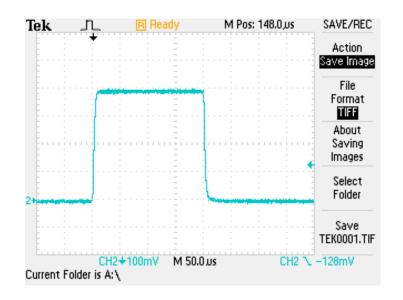

Reference 3000 Booster Current Pulses on 200 m Ω

Here (**Fig. 5**) Gamry Instruments tested a range of current pulses. The smallest current of 2 A did not require the 30k Booster. The larger currents (4 A through 12 A) included the Booster. The curves show no evidence of slew-rate limiting.

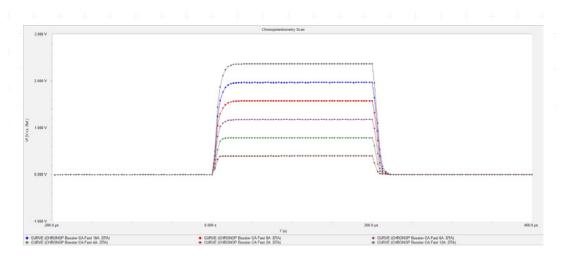


Fig. 5. Data from a range of currents. The lowest current (2 A) used no 30k Booster; the higher currents (4–12 A) added the Booster.

Fig. 6 shows the effect of CA speed on 8 A pulses. The blue curve was recorded with CA speed **fast** and red curve was recorded with CA speed **fastest**.

The red curve has a squarer response but ringing is seen on both the rising and falling edges.

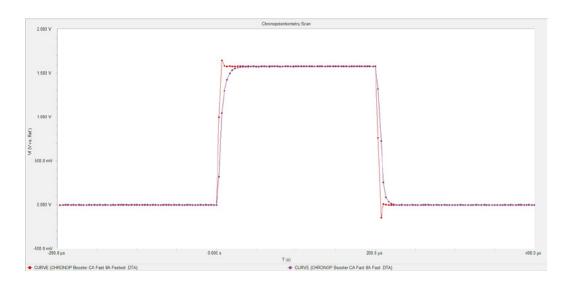


Fig. 6. Data at 8 A, using the CA speed fast (blue) and CA speed fastest (red).



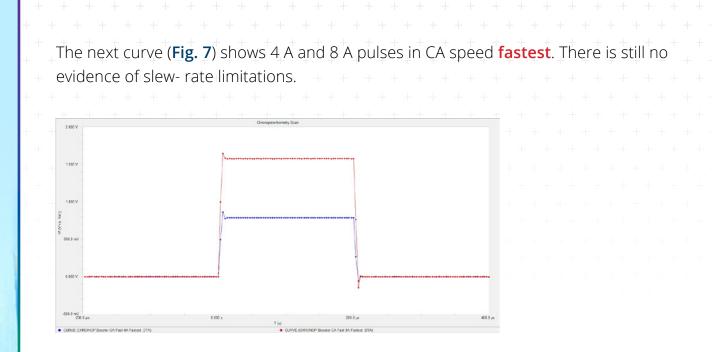


Fig. 7. Data using the CA speed **fastest**, comparing signal responses at 4 A (blue) and 8 A (red).

Results on a 3 m Ω Cell

The connections were switched to the Test side of the Test and Cal Cell. This side is a 3 m Ω , 4 W resistor. At 30 A, its power dissipation is only 2.7 W, so it is safe from misconnection and oscillation.

Fig. 8 shows 2 A pulses on this cell. Such current pulses always are applied with the Reference 3000. Voltage is on the top set of traces and current is on the bottom set. The trace colors correspond to CA speeds: green is **fastest**, red is **fast**, and blue is **norm**.

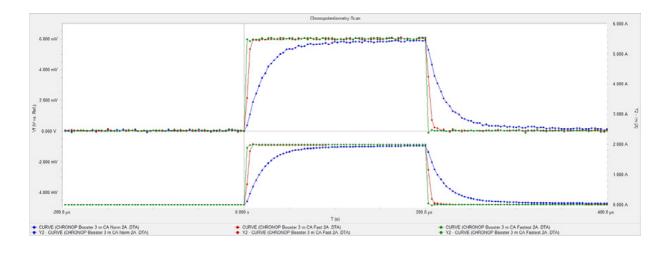


Fig. 8. Data on the Test side of the Test and Cal cell, from 2 A pulses, at CA speeds of **norm** (blue), **fast** (red) and **fastest** (green).

There is a slight indication of ringing in the fastest CA speed.



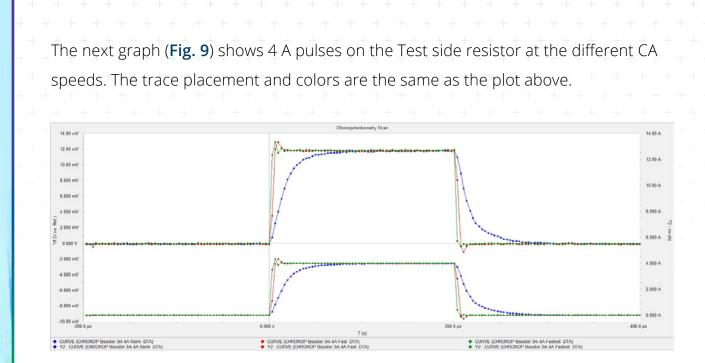


Fig. 9. Data on the Test side of the Test and Cal cell, from 4 A pulses, at CA speeds of **Norm** (blue), **fast** (red) and **fastest** (green). Top: Voltage; bottom: Current.

Both **fastest** and **fast** CA speeds show ringing in **Fig. 9**. The oscilloscope capture in **Fig. 10** shows the **E Monitor** waveform for CA speed **fastest** and a 4 A pulse. Ringing is also visible here.

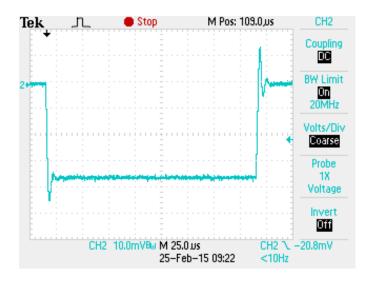


Fig. 10. Oscilloscope data on the Test side of the Test and Cal cell, from 4 A pulses, at the CA speed of **fastest**.



8

50 µs Current Pulses

Fig. 11 shows 50 μ s 2 A and 4 A current pulses into the 200 m Ω Calibrate side of the Test and Calibrate cell. The green trace is a 4 A pulse in CA speed **Fastest**, the red trace is a 4 A pulse with CA speed **fast**, and the blue trace is a 2 A pulse at CA speed **Fast**.

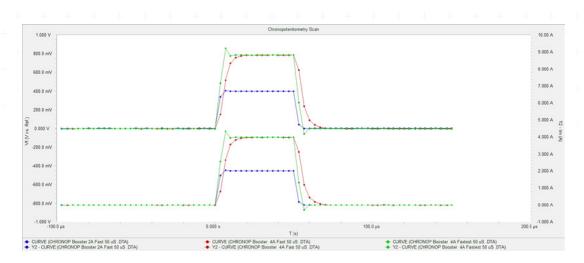


Fig. 11. Data on the Calibrate side of the Test and Cal cell, from 50 µs pulses, at 2 A (blue), 4 A **fast** (red) and 4A **fastest** (green). Top: Voltage; bottom: current.



Summary

Gamry's Reference 3000 and Reference 30k Booster can accurately and reliably generate current pulses as short as the high tens of microseconds.

As to which is the optimal CA speed setting, Gamry Instruments recommends that users run a preliminary test that measures the rise time on a typical cell prior to running a set of experiments.

Galvanostats generally become more stable when cells are capacitive. This effect was not measured in this application note.

Instrument	Cell	τ for Fastest	τ for Fast	τ for Norm
Reference 3000	200 mΩ	τ < 6 μs ringing	τ≈10 µs	τ ≈ 29 μs
	3 mΩ	τ < 6 μs ringing	τ≈7 μs	τ ≈ 23 μs
30k Booster	200 mΩ	τ < 5 μs ringing	τ≈7 μs	τ≈17 μs
	3 mΩ	τ < 2 μs ringing	τ≈6 µs	τ ≈ 22 μs

Approximate rise times τ are shown in the Table below:

Rev. 1 8/17/2017 © Copyright 2017 Gamry Instruments, Inc.



Additional Resources

If you need additional information, please visit our website for our online resources and support.



Technical Support

A compilation of technical support information by hardware and software.



Software Updates

Licensed users can download the latest version of the Gamry Software.



Documentation Downloads Download manuals for hardware and software

products.



Electrochemistry Courses and Training

Information on short courses and electrochemistry training.



Contact Technical Support Get support from the electrochemical experts at Gamry.



Support & Tutorial Videos For all of our tutorial and

support-related videos you can also visit Gamry Instruments YouTube



Application Notes

A series of Application and Technical Notes to assist you in getting the most accurate results.



Literature Database

A database of articles where people have used a Gamry System in their research



For help from our technical staff or questions about our products, please call us at 215-682-9330.





734 Louis Drive • Warminster, PA 18974 • Tel. 215 682-9330 Fax 215 682-9331 • gamry.com